



Transferring DataMate™ to the TI-73 Explorer™

The DataMate software comes already loaded on the CBL 2™. Before you use the CBL 2 for the first time, you must transfer the DataMate software to the TI-73 Explorer.

To transfer DataMate to a TI-73 Explorer, follow these steps:

1. Connect the TI-73 Explorer to the CBL 2 with the I/O init-to-unit link cable.
2. Put the calculator in RECEIVE mode.
 - a. Press **[APPS]**.
 - b. Press **[▶]** to select RECEIVE
 - c. Press **[ENTER]**.
3. On the CBL 2, press TRANSFER. The program or App is transferred and appears in the calculator's list or application list.
4. When the transfer is complete, press **[2nd][QUIT]** on the calculator.

Need Help?

Contact Texas Instruments:

<http://education.ti.com>

ti-cares@ti.com

1-800-TI-CARES